

VALIDIERUNG EINES IMMUNOAFFINITÄTS
LC-MS/MS ASSAYS FÜR DIE SPEZIES-
UND GEWEBESPEZIFISCHE
QUANTIFIZIERUNG VON VERBOTENEN
VERARBEITETEN RUMINANTEN
PROTEINEN IN FUTTERMITTELN


Bundesinstitut für Risikobewertung
Dr. Oliver Pötz
23-Nov-2023





COMPANY

- Founded 2016, fifteen employees, three mass specs
- Novel immunoaffinity mass spectrometry platform for protein quantification
- Industry acceptance: Eight clients from top 20 pharma companies
- Scientific proof: More than 30 peer-reviewed publications
- Critical Path Institute: Predictive Safety Testing Consortium (PSTC) collaboration partner
- Innovation Award Baden Wuerttemberg 2019

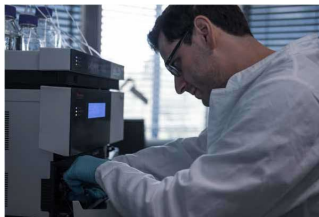

- Challenge – Processed Animal Protein Detection
- Technology – Immunoaffinity LC-MS/MS
- Ring Trial Samples
- Validation
- Summary & Outlook

CHALLENGE – PROCESSED ANIMAL PROTEIN DETECTION

- Difficult samples



- Species identification



- Tissue differentiation



- Sensitive detection & quantification




- Challenge – Processed Animal Protein Detection
- Technology – Immunoaffinity LC-MS/MS
- Ring Trial Samples
- Validation
- Summary & Outlook



SIGNATOPE

TECHNOLOGY – IMMUNOAFFINITY-LC-MS/MS (IA-LC-MS/MS)

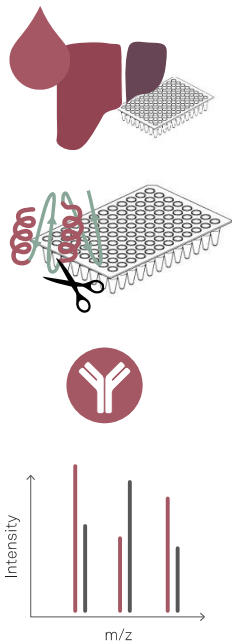

- Sample
- Protein digest using trypsin down to peptides
- Add isotope-labelled peptide standards
- Enrich peptide standards and endogenous peptide derived from protein of interest using antibodies
- Analysis using multiplex targeted nLC-MS/MS

Quantification by ratio of endogenous peptide : internal isotope-labelled standards

SPECIES IDENTIFICATION ASSAY


Heterogeneous
Phase Digestion



Cross-Species
Immunoenrichment




LC-MS/MS
Identification & Quantification

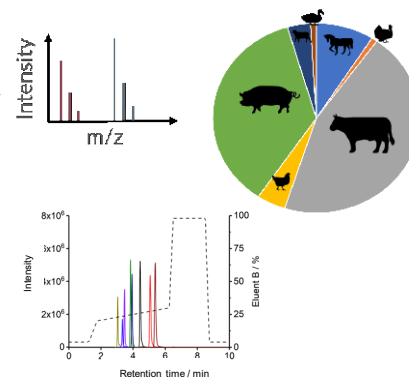

Steinhilber, A. E. *et al. J. Agric. Food Chem.* 2018, 66, 39, 10327-10335

RUMINANT TISSUE DIFFERENTIATION ASSAY

Heterogeneous
Phase Digestion

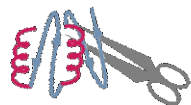



Immunoenrichment
Tissue-Specific Peptides

A2M
SERPINF2
HP252
C9
MYH7
SPP1
MATN1




+ SIS

LC-MS/MS
Identification & Quantification

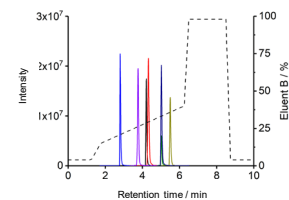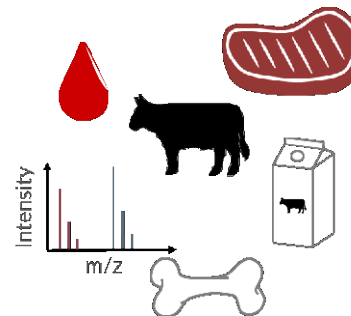


FEED ANALYSIS – A TWO-TIER APPROACH

1. Multi-Species Detection
„which species are present?“

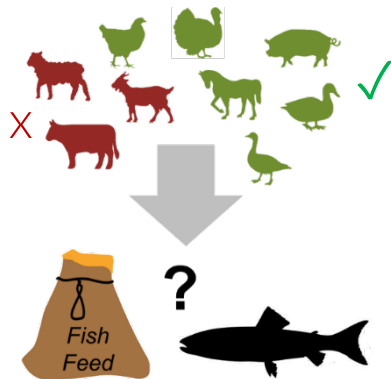

Multiplex 1

2. Tissue-Specific Ruminant Detection
„legal or not?“

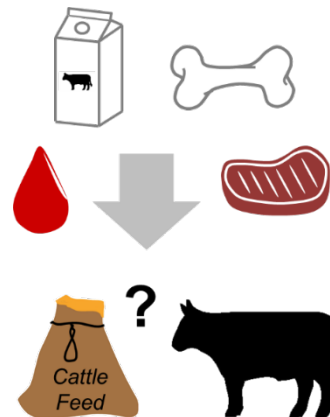

Multiplex 2

- Challenge – Processed Animal Protein Detection
- Technology – Immunoaffinity LC-MS/MS
- Ring Trial Samples
- Validation
- Summary & Outlook











RING TRIAL SAMPLE ANALYSIS

Ring trial samples from former proficiency tests provided by the European Reference Laboratory for Animal Proteins (EURL-AP), Gembloux, Belgium

Sample ID	Description	Additional information about the matrix
Feed 1	0.1 % Ruminant PAP	Pig feed
Feed 2	0.1 % Ruminant PAP	Pig feed
Feed 3	1 % Ruminant Blood	Trout feed, compound feed for trout farming
Feed 4	3 % Bovine Plasma	Trout feed, complete feed for fry
Feed 5	5 % Porcine Blood	Salmon feed
Feed 6	Feed containing hemoglobin meal	Fish feed








SPECIES IDENTIFICATION TIER 1





	Species/ Tissue/ Analyte	Mean peptide amount / fmol						
		Feed 1	Feed 2	Feed 3	Feed 4	Feed 5	Feed 6	Control
Species Identification	Cattle 	4.1 ± 0.1	4.4 ± 0.4	88.0 ± 2.9	956.9 ± 32.3	0	0	0
	Sheep/Goat 	0	0	0	0	0	0	0
	Pig 	0	0	834.3 ± 29.7	0	136.8 ± 4.2	537.0 ± 4.7	0
	Horse 	0	0	0	0	0	0	0
	Turkey 	0	0	0	0	0	0	0
	Chicken 	0	0	0	0	0	0	0
	Duck 	0	0	0	0	0	0	0
	Goose 	0	0	0	0	0	0	0

→ Feed 1-4 are positive for bovine A2M and/or porcine A2M








→ Legal or illegal sources?

TISSUE DIFFERENTIATION TIER 2

	Species/ Tissue/ Analyte	Mean peptide amount / fmol						
		Feed 1	Feed 2	Feed 3	Feed 4	Feed 5	Feed 6	Control
Ruminant tissue Differentiation	SERPINF2 	0	0	2.3 ± 0.1	156.3 ± 7.0	0	0	0
	C9 	0.4 ± 0.1	0.4 ± 0.05	1.2 ± 0.1	50.1 ± 2.5	0	0	0
	HP252 	1.2 ± 0.2	1.4 ± 0.1	10.7 ± 0.3	331.5 ± 10.0	0	0	0
	A2M 	4.5 ± 0.3	4.7 ± 0.5	94.0 ± 4.3	1065.1 ± 165.9	0	0	0
	SPP1 	0	0	0	0	0	0	0
	MYH7 	5.6 ± 0.4	6.7 ± 1.0	0	0	0	0	0
	MATN1 	7.8 ± 1.5	14.4 ± 4.3	0	0	0	0	0

						
	Pig Feed +	Pig Feed +	Trout Feed +	Trout Feed +	Salmon Feed +	Fish Feed +
Samples	0.1 % Bovine PAP	0.1 % Bovine PAP	1 % Bovine Blood + porcine blood (!)	3 % Bovine Plasma	5 % Porcine Blood	containing Hemoglobin Meal


TISSUE DIFFERENTIATION TIER 2

Protein	Sequence	Tissue	Peptide amount / fmol	C.V. / %
A2M	GSGGTAEHPTVEEFVLPK		4.7	11.4
SERPINF2	LPPLSLLK		0.0	-
C9	YTPVEAIEK		0.4	10.9
HP252	FGFDIELFQHAVK		1.4	4.4
SPP1	YPDVAVATWLKPDPSQK		0.0	-
MYH7	MLSSLFANYAGFDPIEK		6.7	14.2
MATN1	AGGIELFAIGVGR		14.4	29.4

→ Feed 2 contains ruminant blood, **meat and bone proteins**

→ no SPP1 detected, hence no milk

→ quantified on a level of 0.1%

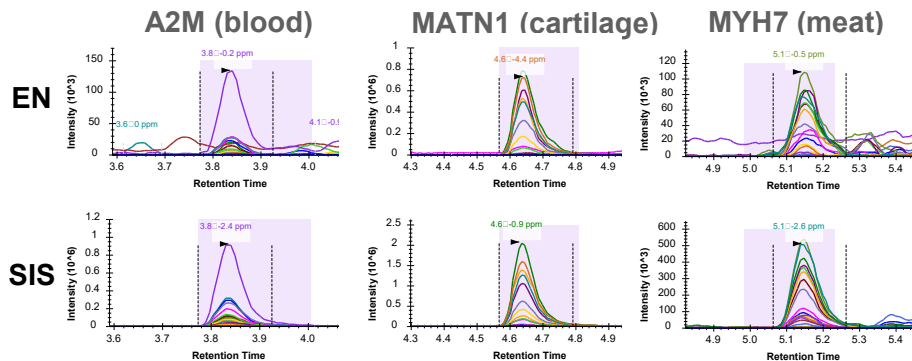



SIGNATOPE

- Challenge – Processed Animal Protein Detection
- Technology – Immunoaffinity LC-MS/MS
- Ring Trial Samples
- **Validation**
- Summary & Outlook









SIGNATOPE

VALIDATION – LOWER LIMIT OF QUANTIFICATION

- Accuracy & Precision – calibration curves & 4 quality controls (0.1, 0.5, 1, 5 % animal protein)
- Limit of Quantification – dilution of calibrators in vegetal feed
- Carry over
- Parallelism – dilution of MBM and SDP
- Selectivity – vegetal feed
- Interference – milk powder
- Analyte stability QCs and calibrators – freeze/thaw, long-term (RT), auto sampler



VALIDATION – LOWER LIMIT OF QUANTIFICATION

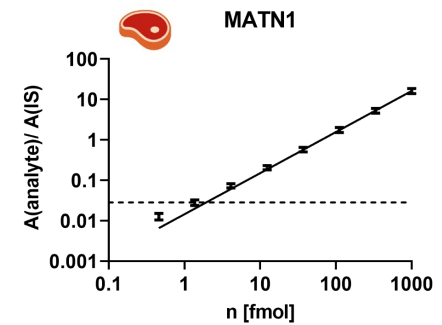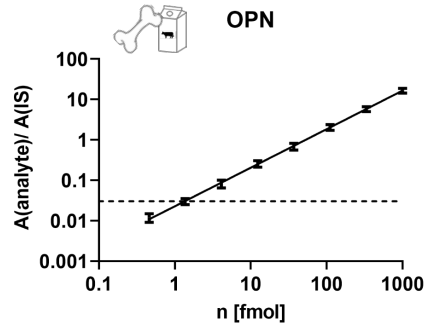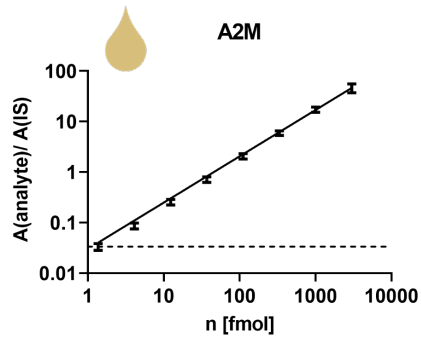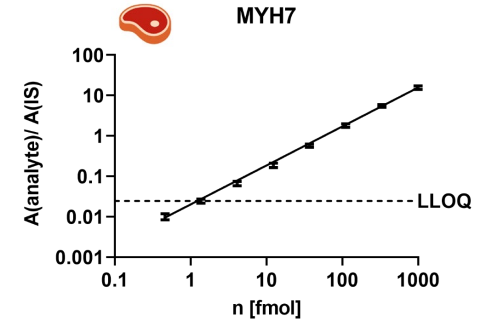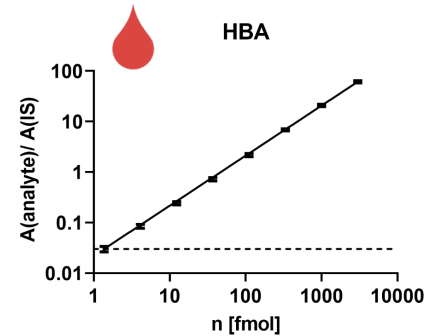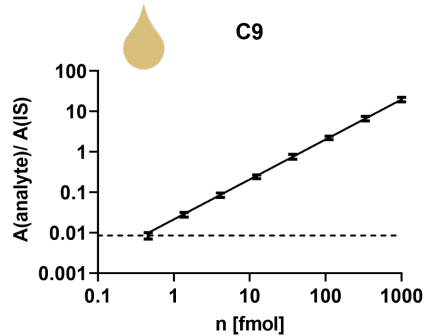

VALIDATION – RESULTS

Parameter	Samples	A2M	C9	MATN1	MYH7	OPN	HBA
LLOQ (ng/g)	Calibrators	15.1	1.82	1.56	20.4	2.67	1.38
ULOQ (ng/g)	Calibrators	33000	3980	3420	14900	1950	3010
Inter precision	Calibrators	passed	passed	passed	passed	passed	passed
Inter accuracy	Calibrators	passed	passed	passed	passed	passed	passed
Inter precision	QC2-QC4*	passed	passed	failed	passed	passed	passed
Inter accuracy	QC2-QC4*	passed	passed	passed	passed	passed	passed
Intra precision	QC2-QC4*	passed	passed	failed	passed	passed	passed
Intra accuracy	QC2-QC4*	passed	passed	passed	passed	passed	passed
Reproducibility	contrieved samples (one digest, IP n=2)	passed	passed	passed	passed	passed	passed
	contrieved samples (digest, n=2)	passed	passed	failed	passed	failed	failed

VALIDATION – RESULTS

Parameter	Samples	A2M	C9	MATN1	MYH7	OPN	HBA
Selectivity	spiked samples / unspiked samples	passed	failed	failed	passed	failed	passed
Parallelism	commercial/artificial samples, serially diluted	passed	passed	failed	passed	passed	passed
Carryover	BS1-BS8	>S2	>S2	>S4	>S3	passed	passed
Interference	spiked samples, spiked milk powder	failed*	failed*	failed*	Up to 60 %	failed*	Up to 10%
Freeze-thaw stability	Proteolyzed sample, 1x-3x	failed	3x	3x	3x	3x	3x
	calibrators, 1x-3x	passed	passed	passed	passed	passed	passed



VALIDATION – SUMMARY

- **Accuracy & Precision:** 0.1% animal protein QC1 failed the quantitative acceptance criteria but not the qualitative acceptance criteria.
- **Limit of quantification:** Method sensitive to detect PAP down to 0.1%.
- **Reproducibility:** Technical measurement passed acceptance criteria, but 3 out of 6 assays failed most probably due to sample homogeneity.
- **Carry over:** Samples with content > S3 must be re-run with additional blank run
- **Parallelism:** Passed, but not MATN assay.
- **Selectivity:** A2M, MHY7 and HBA assays met criteria.
- **Interference:** Only MHY7 and HBA assays met criteria, milk powder contains OPN, C9, A2M & MATN.
- **Analyte, QCs and calibrators:** Long-term stability confirmed for up to 30 months.
- **Reagent stability:** Confirmed for up to 30 months.



SIGNATOPE

- Challenge – Processed Animal Protein Detection
- Technology – Immunoaffinity LC-MS/MS
- Ring Trial Samples
- Validation
- Summary & Outlook

SUMMARY & OUTLOOK

- Method sensitive to detect PAP down to 0.1% & to quantify down to 0.5%
- Sampling process requires optimization, sample size needs to be determined to achieve statistical robust result
- Interference: MHY7 tolerated up to 60% milk powder
- Reagent stability: Confirmed for up to 30 months
- Ring trial in progress
 - Dr. Uta Herfurth German Federal Institute of Risk Assessment, Germany
 - Prof. Rene Zahedi, University of Manitoba, Canada
 - Dr. Ikram Belgit, Institute of Marine Research, Norway
 - Dr. Marie Lecrenier, Wallon Agriculture Research Centre, Belgium
 - Prof. Jens Brockmeyer, Universität Stuttgart, Germany

TEAM & FUNDING



Oliver Pötz
Hannes Planatscher
Thomas Joos
Isabel Bek
Katharina Bendel
Sonja Burger
Helen Hammer
Lena Heumesser
Hanna Hitzelberger
Robin Kretz
Tobias Meisinger
Wael Naboulsi
Anja Tausch
Amelie Vogt
Cornelia Sommersdorf

